

Generating a Devyser LB-192 Plate sample sheet with Illumina LRM

Generate a sample sheet for each Devyser run by using the Illumina Local Run Manager (LRM) software according to the current software guide³. The LRM can be installed on the Illumina® MiSeq™, MiniSeq™ or iSeq™ 100 computer.

Before proceeding with these instructions, ensure that the LRM 'GenerateFASTQ' Analysis Module is installed.

These instructions are applicable for LRM v2 and later^{1,2,3}. Please contact techsupport@devyser.com if the software version is incompatible.

1. LRM File Import

Download the zip file 'Index Plate LB-192 LRM files vYYYY-MM-DD' from the Devyser website and unzip the files.

- A. Open LRM and select 'Library Prep Kits' from the TOOLS drop-down menu
- B. Select 'Add Library Prep Kit'
- C. Import the un-zipped .tsv file (Devyser 2x76 LB-192, Devyser 2x151 LB-192 or Devyser 1x301 LB-192 file) and ensure that it is present in the Library Prep Kits list

2. Sample sheet generation

- A. From the RUN DASHBOARD, select 'Create Run' and then select 'GenerateFASTQ'
- B. Enter a 'Run Name' and 'Run Description' of your choosing
- C. Select the previously imported Library Prep Kit from the Run Settings subheader
- D. Scroll down and select the '+' symbol besides 'Rows' to add the desired number of rows/samples
- E. Enter the sample names in the Sample ID column
Note: When using a MiSeq, select the plate-location of each sample in the INDEX WELL column. Go to step H
- F. In the INDEX 1 (I7) column, choose the correct index (LB001-LB192) in the drop-down menu (see table 1)
- G. In the INDEX 2 (I5) column, choose the correct index (LB001-LB192) in the drop-down menu (see table 1)
- H. Click on 'Save Run', or 'Export Sample Sheet' to be imported at a later date

3. Index Description

The Illumina double indexes introduced during PCR2 are listed in table 1 and 2.

Table 1: Illumina double indexes used in Index plate LB-1

	1	2	3	4	5	6	7	8	9	10	11	12
A	LB001	LB009	LB017	LB025	LB033	LB041	LB049	LB057	LB065	LB073	LB081	LB089
B	LB002	LB010	LB018	LB026	LB034	LB042	LB050	LB058	LB066	LB074	LB082	LB090
C	LB003	LB011	LB019	LB027	LB035	LB043	LB051	LB059	LB067	LB075	LB083	LB091
D	LB004	LB012	LB020	LB028	LB036	LB044	LB052	LB060	LB068	LB076	LB084	LB092
E	LB005	LB013	LB021	LB029	LB037	LB045	LB053	LB061	LB069	LB077	LB085	LB093
F	LB006	LB014	LB022	LB030	LB038	LB046	LB054	LB062	LB070	LB078	LB086	LB094
G	LB007	LB015	LB023	LB031	LB039	LB047	LB055	LB063	LB071	LB079	LB087	LB095
H	LB008	LB016	LB024	LB032	LB040	LB048	LB056	LB064	LB072	LB080	LB088	LB096

Table 2: Illumina double indexes used in Index plate LB-2

	1	2	3	4	5	6	7	8	9	10	11	12
A	LB097	LB105	LB113	LB121	LB129	LB137	LB145	LB153	LB161	LB169	LB177	LB185
B	LB098	LB106	LB114	LB122	LB130	LB138	LB146	LB154	LB162	LB170	LB178	LB186
C	LB099	LB107	LB115	LB123	LB131	LB139	LB147	LB155	LB163	LB171	LB179	LB187
D	LB100	LB108	LB116	LB124	LB132	LB140	LB148	LB156	LB164	LB172	LB180	LB188
E	LB101	LB109	LB117	LB125	LB133	LB141	LB149	LB157	LB165	LB173	LB181	LB189
F	LB102	LB110	LB118	LB126	LB134	LB142	LB150	LB158	LB166	LB174	LB182	LB190
G	LB103	LB111	LB119	LB127	LB135	LB143	LB151	LB159	LB167	LB175	LB183	LB191
H	LB104	LB112	LB120	LB128	LB136	LB144	LB152	LB160	LB168	LB176	LB184	LB192

4. References

- ¹Local Run Manager v2 Software Guide (Illumina document # 1000000002702)
- ²Local Run Manager v3 Software Guide (Illumina document # 1000000111492)
- ³Local Run Manager v4 Software Guide (Illumina document # 200036672 v00)

5. Revision History

Version 2026-06-10

New.